



Illumina Proactive Service connectivity guide

Easily connect your NovaSeq™ 6000 Sequencing System

To reduce unplanned downtime, minimize run failures and lost samples, plus resolve issues faster if and when they arise, simply follow the steps below.

Follow these steps to get connected

- From the pull-down menu on the **Home** Screen, select **Settings**
- Under the **Mode Selection** tab, ensure the box next to **Send instrument performance data to Illumina** is selected
- Click **Save**

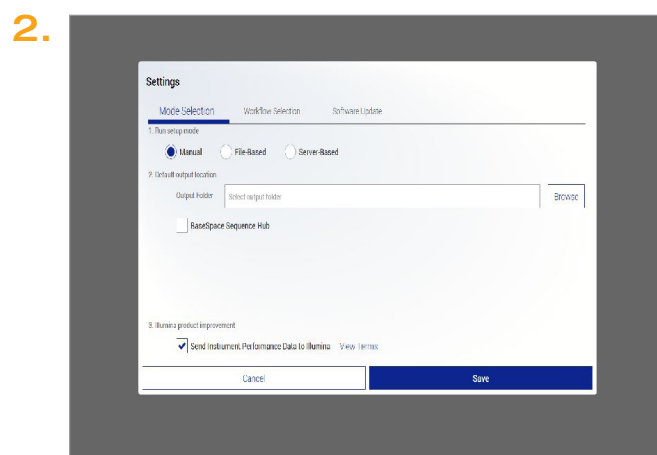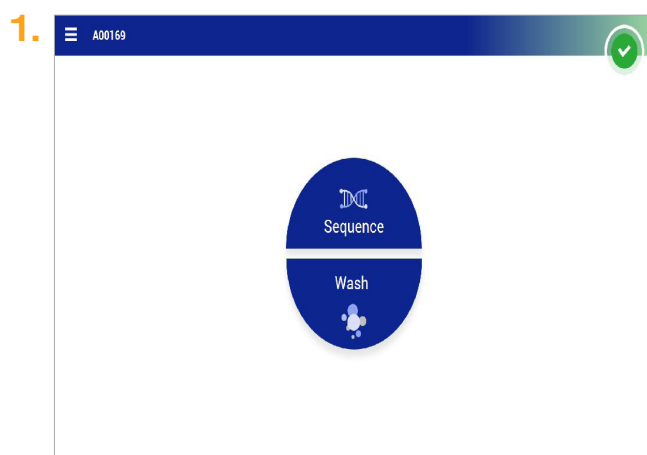

For more information, contact your local field applications scientist or field service engineer.

www.illumina.com

For Research Use Only. Not for use in diagnostic procedures.